

POLYTRON® PT 3100 D

Stand-disperser (High-End Line)

FLEXIBILE-USE DISPERSER

With digital controller, numerous applications « Working volume 0.1 to 10,000 ml »

R

POLYTRON PT 3100 D





POLYTRON® - Efficient, Powerful, Digital Controls

Applications with the highest demands for reproducibility and scaling up can be efficiently performed with this device. The fully digitized control and regulation system is very easy to operate. The illuminated LCD display shows all the parameters, which can be changed with the membrane keypad.

TECHNICAL INFORMATION

Processing volume (water) 0.1 to 10,000 ml

Tip speed Up to max. 32 m/s

Applicable aggregates ø 5 to 36 mm

Speed range 500 to 30,000 rpm

Drive coupling Aggregates with F coupling

Noise level (drive without aggregate) 66 dB(A) at 25,000 rpm 73 dB(A) at 30,000 rpm

Motor Universal with carbon brushes

Drive power 1200W

Connection voltage 110 - 230 V ± 10%, 50Hz / 60Hz

Relative humidity (max.) 80% during storage 80% during operation

Operating temperature 0 – 40 °C

PC interfaces RS-232 (D SUB 9) USB socket B

Protection class acc. to DIN IP 20

Dimensions (L x W x H) Drive: 240 x 130 x 265 mm

Weight Drive: 3.9 kg

EMC standards IEC/EN 61000-6-2 / EN 61000-6-3

Safety standard IEC/EN 61010-2-51



THE EXPERT FOR YOUR LABORATORY

For more than 60 years, we have been perfecting the rotor/ stator technology, working in close cooperation with users and universities. When using our technology, your test samples will be dispersed into homogeneous, chemical systems. Depending on the type of samples you will get homogeneous emulsions, dispersions or suspensions. This is the best foundation for all subsequent analyses or formulations.

ERGONOMICS

- Easy operation with an on-off switch and membrane keypad
- Digital speed control for long equipment life
- Protection against overheating, overload and blocking
- Speed stability even with changes in viscosity
- Easy to adjusted motor height (depending on the stand)
- Aggregates guickly switched with Click & Go coupling
- Easy aggregate clean–up with EasyClean Design

SCALE-UP



You can use test results to scale-up to our pilot plant and industrial POLYTRON[®] machines. The volume range for on batch units is seamlessly available up to 10,000 liters.

RESULTS ARE CONVINCING

SMALLEST SIZES OF PARTICLES/DROPLETS because precise analysis require homogeneous samples

EFFICIENT SAMPLE PREPARATION because fast dispersion saves time

STANDARD OR CUSTOM-MADE AGGREGATES perfected for and in cooperation with the users

CLEANING AND DISASSEMBLY EasyClean Design makes it very simple


QUICK ENGAGEMENT OF THE AGGREGATES by using Click & Go

BEST MATERIALS AVAILABLE because without them, we could not call it Swiss Made

DURABILITY is a Swiss virtue



Users / Application ranges

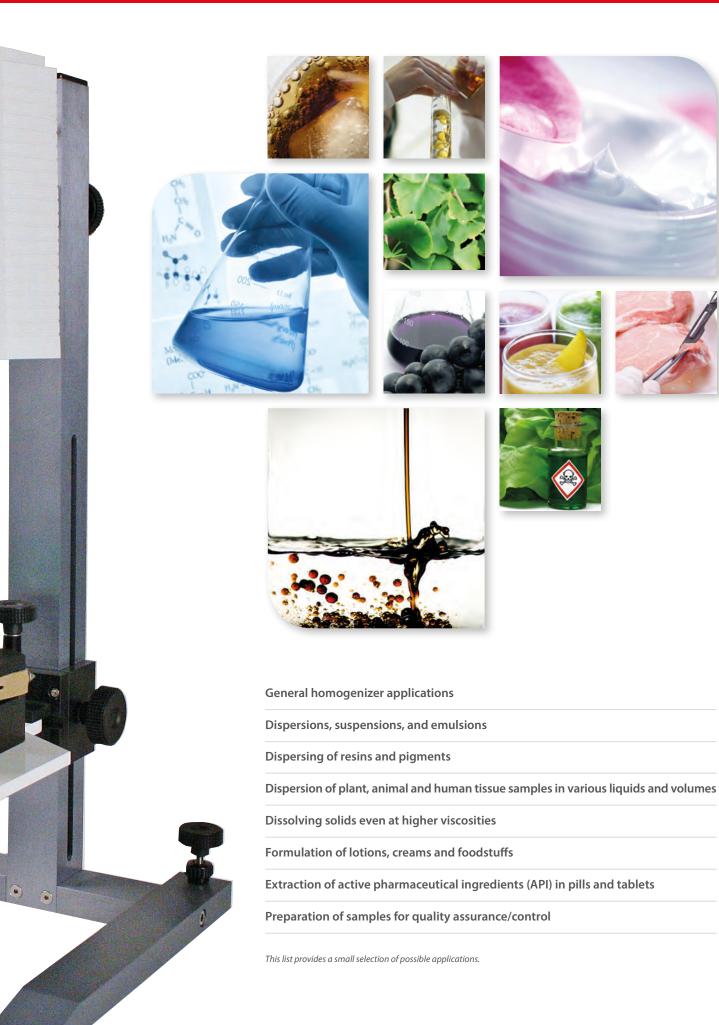

Kinematic dispersion using the rotor/stator systemtem

Who is the founder of it?

The roots of today's KINEMATICA began in the former chemical/physical research institute of Prof. P. Willems in Lucerne Switzerland, the inventor of the modern rotor/stator dispersion technology.

THE PRINCIPLE:

The spinning rotor generates a vacuum. The sample is drawn in and discharged to the outside through the stator's slots. Between the rotor/stator (shear gap), the product is subject to high deceleration-tangential and radial acceleration forces. The individual particles/droplets are torn apart and thus reduced in size; they are reduced several micrometers in size through additional cutting and crashing effects.

THE RESULT:

Microscopic dispersions, emulsions, suspensions, and foams are generated. Droplets, particles, and gas bubbles are reduced to several micrometers or smaller. Small and large substances are dispersed more economically, faster, and better than with any other system.

Your samples are efficiently «POLYTRONIZED».

Aggregate Technology

Unrivaled!

The high efficiency and the uniqueness of our POLYTRON[®] dispersing aggregate can only be achieved through superior manufacturing precision and the sophisticated KINEMATICA design method. This is 100% Swiss Made.

STANDARD AGGREGATES (EC DESIGN, WITH TWO OR MORE SPROCKETS)

This is all you need for popular applications in the preparation of samples and formulation development. The cutting teeth embedded in the head of the tool optimize the pre-crushing of the sample. An enclosed ring design provides optimum protection for those teeth and prevents them from bending.

SPECIAL AGGREGATS (W, M, B, DS, Z DESIGN)

Specializations in the chemical, pharmaceutical, cosmetic, and food industry, as well as the life sciences, makes new designs inevitable. These aggregates are based on state-of-the-art technology and only available from KINEMATICA.

The highly specialized tool is subject to further development that involves close cooperation with our customers and universities. Thus, it is continuously perfected for specific applications.

MATERIALS

Whether you need a tool with standard or special geometry, all aggregates are manufactured from high-alloy stainless steel 316L, electrolytically polished and made in Switzerland.

BEARINGS


Depending on the diameter, our aggregates are made with high-quality plastic, special carbide or bronze/carbide bearings. They have excellent mechanical properties and chemical resistance and have FDA approvals. Super high-speed ball bearings are used in the couplings to provide stability at speeds of up to 30,000 rpm.

CLEANING AND DISASSEMBLY

EasyClean: Our technology ensures the simplest removal and fast cleaning, either by hand, dishwasher, or autoclave. Disassembling the precisely fitting components is simple and fast.

LOCKING SYSTEM

All our POLYTRON® aggregates are fitted with a high-precision quick-release system: The Click & Go

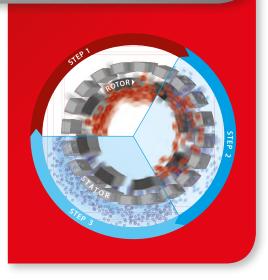

Dispersing aggregates for PT 3100 D

Your applications, our response

KINEMATICA aggregates solve all problems you are faced with every day in the laboratory. Using our Click & Go quick release makes the changing of aggregates very easy.



G «GAS TIGHT»

EC «STANDARD DISPERSING AGGREGATE»

Design



The universal geometry uses two rows of teeth, is suitable for a large spectrum of laboratory applications, and it is reliable and efficient

- Sample preparation for medical diagnostic procedures as well as DNA analysis - Extractions of systems/substances
- from vegetable-derived samples - Disruption of animal and human cells

- Deagglomeration of united cell

structures

- Selection and applications
- All standard dispersing taskes - Grinding of animal and human tissue samples in a variety of fluids - Intensive mixing and volumes
- Sample preparation for the extraction and dissolution of organic materials

Sales No.:	Description	ø/length mm	Volume ml	max. Tip speed m/s
11030297	PT-DA 05/2EC-F078	5 / 78	0.1 - 5	5
11030298	PT-DA 07/2EC-F101	7 / 101	0.3 - 10	8
11030299	PT-DA 12/2EC-F154	12/154	3 - 250	14
11030302	PT-DA 20/2EC-F193	20 / 193	20 - 2'000	22
11030305	PT-DA 30/2EC-F250	30 / 250	100 - 4'000	30
11030311	PT-DA 30/4EC-F250	30 / 250	100 - 4'000	30
11030313	PT-DA 36/2EC-F250	36 / 250	200 - 4′500	34
11030317	PT-DA 36/4EC-F250	36 / 250	200 - 4′500	34

	M «OUTSI	DE CUTTING BLA	DES»			
All and		Design Outside blades pre-c that are larger than t diameter. Subsequer be dispersed inside t	rush samples he rotors ntly, they can			
	Sales No.:	Description	ø/length mm	n Volume ml	max. Tip speed m/s	5

Sules No.	Description	ø/iengarinin	volume im	max. np speed m/s
11030300	PT-DA 12/2MEC-F154	12 / 154	3 - 250	14
11030303	PT-DA 20/2MEC-F193	20 / 193	20 - 2′000	22
11030307	PT-DA 30/2MEC-F250	30 / 250	100 - 4'000	30
11030308	PT-DA 30/2MG-F273*	30 / 273	100 - 4'000	30
11030315	PT-DA 36/2MEC-F250	36 / 250	200 - 4'500	34

* M-design aggregate in gas-tight model

W «**W-DESIGN**»



Design The W design prevents fibrous, stringy or solid samples from clogging the aggregate

Selection and applications - Dispersing fibrous, stringy or solid samples (e.g., pieces of meat) - All standard dispersing applications similar to EC aggregate

Sales No.:	Description	ø/length mm	Volume ml	max. Tip speed m/s
11030301	PT-DA 12/2WEC-F154	12/154	3 - 250	14
11030304	PT-DA 20/2WEC-F193	20/193	20 - 2'000	22
11030294	PT-DA 20/2G-F210*	20/210	20 - 2'000	30
11030309	PT-DA 30/2WEC-F250	30 / 250	100 - 4'000	30
11030345	PT-DA 30/2WMEC-F250**	30 / 250	100 - 4'000	30
11030319	PT-DA 36/2WEC-F250	36 / 250	200 - 4'500	34
11030346	PT-DA 36/2WMEC-F250**	36 / 250	100 - 4'500	34

W-design aggregate in gas-tight model ** W-Design aggregate with knifes

Design Integrated	mechanical seals	Selection and ap – Under pressure – Preventing air f added to emuls – Corrosive acid r	or vacuum rom being sions	S.
Sales No.:	Description	ø/length mm	Volume ml	max. Tip speed m/s
11030295	PT-DA 20/2G-F210	12/210	20 - 2'000	22
11030306	PT-DA 30/2G-F273	30/273	100 - 4'000	30
11030312	PT-DA 30/4G-F273	30/273	100 - 4'000	30
11030316	PT-DA 36/2G-F273	36/273	200 - 4'500	34
11030318	PT-DA 36/4G-F273	36/273	200 - 4′500	34

				Z « Z-DESIGN »	
integrated	o of hard samples, with V-notch technology. integrated holes for	Selection and ap – Hard and brittle – Sweet corn and – Breaking, open of pills – Breaking, hard g	material grain and dispersing	ALL	-
Sales No.:	Description	ø/length mm	Volume ml	max. Tip speed m/s	
11030310	PT-DA 30/2ZEC-F250	30 / 250	100 - 4'000	29	

B «BIOTRONA® HIGH-TURBULENCE MIXER»						
Design High-turbulence mixing head with the least possible shear forces and energy inputs		Selection and applications – Quick removal and suspension of solid particles; also suitable for high viscosity applications		10.3		
Sales No.:	Description	ø/length mm	Volume ml	max. Tip speed m/s		
11030296	PT-DA 20/BEC-F193	20 / 193	100 - 2'000	21		
11030314	PT-DA 36/BEC-F250	36 / 250	500 - 10'000	34		

	DS <	«DISSOLVER»
Design A processing head with teeth. T product is radially accelerated	Selection and applications – Fast mixing and homogenizing of solids in liquid – Processing of highly viscous products	-
Sales No.: Description	ø/length mm Volume ml	max. Tip speed m/s
11030344 PT-DA 30/DS-F25	30 / 250 500 - 10'000	30

	TIN « TITANIUM NITRATED» - UPGRADE
Design Rotor and/or stator with up to three rows of teeth with a titanium nitrated (TiN) coating	Selection and applications - All dispersing applications with abrasive samples
Sales No.: Description	

11035050	TIN rotor for the selected dispersing aggregate, ø 12-36 mm
11035150	TiN stator for the selected dispersing aggregate, ø 12-36 mm

Dispersing vessels

Choosing the right dispersing aggregate is crucial for good results. The right vessel also helps contribute to achieving the desired results.

THE PROBLEM

For normal cylindrical beakers, the flow creates a vortex. The mixing and dispersing efficiency decreases considerably since the product is no longer being optimally processed through the dispersing generator head. This lengthens processing times, requiring more energy input. Depending on the shape of the vortex, this may also lead to additional air being introduced.

THE SOLUTION

Our specially designed GS jars have a cloverleaf-like shape. This effectively interrupts the flow in the radial direction and generates additional counteracting forces. These high turbulences maximize the mixing and dispersion efficiency, thus minimizing processing time and requiring less energy.

THE SELECTION

Our GS-glasses are available in chemical- and temperature-resistant borosilicate glass or in stainless steel, with working volumes from a few milliliters to several liters, with or without lid and with or without sealed feedthrough.



GS vessel principle (viewed from above)






PC control software


CONTROL PANEL

The PT 3100 D is controlled digitally via a keypad. All important operating parameters can be read on the 2-line LCD display

CE OFF	ON / OFF switch
POLYTRON	LCD display standard, timer and automatic modes change modes change display settings cursor left/right confirm input values start motor/release control lower/increase value/speed stop motor

The free-of-charge software «User Interface PT 3100 D» was developed for the operation and control of the laboratory dispersing system POLYTRON® PT 3100 D. A USB or RS232 interface provides the communication to the computer. The data transmission protocol complies with the Modbus specifications.

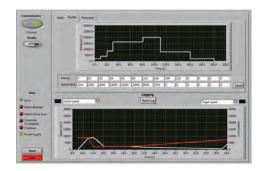

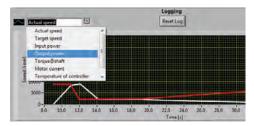

OPERATING MODE

- STANDARD MODE: Operation by manual control
- TIMER MODE: Automatic shutdown by recording speed and time
- AUTO MODE: Complete control through the included KINEMATICA software

KINEMATICA SOFTWARE

- Set the speed via PC
- Store speed/time profile to reproduce formulas or processes
- Adjustable temperature limits as stop criterion, for example, to prevent product overheating (with Pt100 temperature probe)
- Graphical representation of parameters with time diagram
- Continuous recording of all parameters listed above in MS Excel format
- PC software and cable included

ADDITIONAL FEATURES

 Memory function: remembers the last speed setting and the time when the device shut off

 displays the product temperature by means of the optional Pt100 temperature probe

PARAMETERS RETRIEVABLE VIA LCD DISPLAY

- Actual speed [AS, rpm]
- Target speed [TS, rpm]
- Power in [Pin, W]
- Power out [Pout W]
- Moment of force [M Ncm]
- Motor current [I, A]
- Supply voltage [U, V]
- Product temperature (with optional Pt100 temperature probe) [TP, $^{\circ}$ C]
- Internal temperature controller [Tc, °C]
- Operating hours meter [Oh, h]
- Firmware

SYSTEM REQUIREMENTS:

Operating system:	Windows [®] XP / 7 / 8
Required interfaces:	1 x USB or 1 x RS-232





Comfort at the highest level

Modern laboratories want equipment that is easy to handle. This rugged electric tripod offers a variety of benefits.

REQUIREMENTS

The processing of large sample volumes or viscous products requires larger, higher torque drive units and larger aggregates which are quite heavy. These factors were the reasons for making this stand available to our customers.

Now, what has long been a part of the standard equipment in the pilot plants, is now available for laboratory use.

PRINCIPLE

- Powerful motor mounted onto the telescopic lifting arm
- Durable, robust construction
- High stability due to the massive aluminum base plate

RESULT

- Easy height adjustment by means of two buttons
- Anti-slip mat included
- Available with optional vessel holder





OTHER STANDS AND ORDER NUMBERS

Figure	1	2	3	3	3	3
Sales no.:	11040025	11040032	11040033	11040036	11040037	11040034
Designation	ST-F 10/600	ST-F 13/820	ST-P 14/590+500, 230 V (EU-plug)	ST-P 14/590+500, 230 V (CH-plug)	ST-P 14/590+500, 230 V (UK-plug)	ST-P 14/590+500, 100-120 V
Description	H-foot stand made of anodized aluminum and the height adjustable, swiveling PTFE-coated work table with vessel holder. Weight: 5.5 kg	H-foot "Economy" stand with a column, two cross joints, safety ring and vessel holder. If you want to use this stand with the PT 3100 D, you will need the optional adapter bracket 2	Electric plate stand with lifting telescopic bar for more ease of use. Incl. up/down button and mat to prevent slipping.			
			The electronics are desi EN 60335-1 and UL 962		tures of +5℃ to +40℃ and	were tested per
		Weight: 10.5 kg	Weight: 11 kg			
	1					
					STA	AND ACCESSOR





		STAND ACCESSORIES
Sales no.:	Designation	Description
11045013	Adapter Bracket 2	For the "Economy" stand ST-F 13/820
11036005	Vessel holder	Vessel holder with guide for ST-P 14/590+500

Order information and accessories

Depending on your specific requirements, it is possible to expand the PT 3100D drive with our functional accessories. Our expert personnel will be happy to assist you and answer all the questions you might have.

ADAPTER

Adapters for use with dispersing aggregate with A coupling

Sales No.:	Designation
11035001	Adapter PT-DA / PTA (F/A)

Description

This adapter connects older A-coupling dispersing aggregates such as PT-K / PT-G with this PT 3100 D motor (with F-coupling).



	DRIVES / COMPLETE SYSTEMS
Sales No.:	Description
	Motor
11010120	POLYTRON® PT 3100 D, 230 V / CH-Power cord, PC software, RS-232/USB connection cable
11010121	POLYTRON® PT 3100 D, 230 V / EU-Power cord, PC software, RS-232/USB connection cable
11010122	POLYTRON® PT 3100 D, 230 V / UK-Power cord, PC software, RS-232/USB connection cable
11010123	POLYTRON® PT 3100 D, 100-120 V / Power cord, PC software, RS-232/USB connection cable
	Motor is sold with stand and accessories
11090090	PT 3100 D, 230 V, CH-Power cord, stand ST-F10/600, PC software, RS-232/USB connection cable
11090091	PT 3100 D, 230 V, EU-Power cord, stand ST-F10/600, PC software, RS-232/USB connection cable
11090092	PT 3100 D, 230 V, UK-Power cord, stand ST-F10/600, PC software, RS-232/USB connection cable
11090093	PT 3100 D, 100-120 V / Power cord, stand ST-F10/600, PC software, RS-232/USB connection cable



KINEMATICA solutions for your homogenizing, dispersing, emulsifying, suspending, crushing, wet milling tasks.

Our staff will be happy to advise you. We will be glad to you additional documentation or present the POLYTRON[®] systems to you in person.

YOUR APPLICATION IS OUR FOCUS!

KINEMATICA is a leading manufacturer of dispersing and mixing technology for standard and customized applications in the lab, pilot plant and production areas of pharmaceutical, chemical, food, cosmetic and biotech or life science companies worldwide.

Our POLYTRON[®] batch and MEGATRON[®] In-line Homogenizers are suitable for many applications:

- Dispersing of-non soluble liquids or solids into liquids to finest emulsions/ multiple emulsions or suspensions
- Induction and dispersion of powders into liquids
- Foaming by gas induction into liquids
- Disintegration of tissue samples for preparation in further analysis
- Dispersing of various samples for quality control

We also deliver POLYMIX[®] Micro Dry Grinding Mills and a variety of POLYMIX[®] and MICROTRON[®] Overhead Stirrers and Mixers.

Whatever your application, we are confident that our team of specialists, with over 60 years of experience, will provide the best solution for you.

YOUR SATISFACTION IS OUR GOAL!







Batch dispersing/homogenizing units



High turbulent mixing



Inline dispersing/homogenizing units



Dispersing/homogenizing Reactors



Stirrers and dry Mills



Mixer/Blender



Viscometers









Headquarters KINEMATICA AG Dispersing and Mixing Technology

Luzernerstrasse 147a 6014 Luzern Switzerland

 Tel.
 +41 41 259 65 65

 Fax
 +41 41 259 65 75

 E-Mail
 info@kinematica.ch

Subsidiaries KINEMATICA, INC. Dispersing and Mixing Technology

155 Keyland Court Bohemia, NY 11716 USA

 Tel.
 +1 631 750 66 53

 Fax
 +1 631 750 66 57

 E-Mail
 info@kinematica-inc.com

KINEMATICA AG Dispersing and Mixing Technology

Office Eastern Europe Prosp. Akad. Palladina 44 UA – Kyiv 03680

Tel./Fax+38 044 422 61 27E-Mailinfo@kinematica.com.uawebwww.kinematica.com.ua

KINEMATICA GMBH

Dispersing and Mixing Technology

Münstertäler Str. 12 79427 Eschbach Germany

E-

el.	+49 7634 504 800 0
х	+49 7634 504 800 22
Mail	info@kinematica.de