



KINEMATICA

Homogenizing perfected.

POLYTRON®



POLYTRON® SYSTEM PT 10-35 GT

Operating instructions

**Garantie
Warranty
Upgrade**



Jetzt registrieren / Register now:
kinematica.ch/garantie-upgrade
kinematica.ch/warranty-upgrade

OPERATING INSTRUCTIONS			
System	POLYTRON®		
Type	PT 10-35 GT		
Issue	3 / 01.07.2018		
Page	2	of	32

TABLE OF CONTENT:

1 INTRODUCTION	3
1.1 OPERATING INSTRUCTIONS.....	3
1.2 ORGANISATIONAL MATTERS	5
1.3 WARNING NOTICES.....	6
2 SAFETY.....	7
2.1 SUMMARY	7
2.2 SAFETY CONCEPT	7
2.3 RESIDUAL DANGERS.....	10
2.4 WARNINGS.....	11
3 DESCRIPTION OF THE EQUIPMENT	13
3.1 OVERVIEW	13
3.2 DRIVE UNIT.....	14
3.3 STANDS ST-P10/600 & ST-P20/600	16
3.4 POLYTRON® DISPERSING-AGGREGATES	17
3.5 TECHNICAL SPECIFICATIONS	18
4 INSTALLATION	19
4.1 UNPACK.....	19
4.2 ASSEMBLING OF THE DRIVE.....	20
4.3 ASSEMBLING OF THE STAND WITH ACCESSORIES	20
4.4 STARTING UP	21
5 MAINTENANCE	23
5.1 DRIVE	23
5.2 CLEANING AND MAINTENANCE OF THE AGGREGATES.....	24
6 TROUBLE SHOOTING	26
7 ACCESSORIES	27
8 WARRANTY	28
9 DISPOSAL.....	29
APPENDIX A. DIMENSIONAL DRAWINGS	30
APPENDIX B. BILL OF MATERIALS.....	32

OPERATING INSTRUCTIONS			
System	POLYTRON®		
Type	PT 10-35 GT		
Issue	3 / 01.07.2018		
Page	3	of	32

1 INTRODUCTION

This chapter gives information on the the structure of this document. It will assist you in making use of it and show how to find the required information quickly.

1.1 OPERATING INSTRUCTIONS

PLEASE READ THESE OPERATING INSTRUCTIONS BEFORE SWITCHING ON OR OPERATING THE EQUIPMENT. THEY DESCRIBE THE USE OF THE POLYTRON® PT 10-35 GT, ITS INSTALLATION AND MAINTENANCE AND THE APPROPRIATE REPLACEMENT PARTS AND ACCESSORIES.

THEY WILL HELP YOU AVOID ERRONEOUS USE AND SUBSEQUENT DAMAGE. ALTHOUGH POLYTRON® UNITS ARE DESIGNED FOR EASE OF SERVICE, THIS DOES NOT RELEASE YOU FROM THE OBLIGATION TO INSPECT YOUR EQUIPMENT CAREFULLY AND TO CLEAN IT THOROUGHLY.

KINEMATICA AG is a specialist manufacturer of machines and equipment for dispersion and mixing technology.

An important objective of these operating instructions is to fully inform you, the user, about the correct and safe use of our equipment.

In order to achieve this, it is essential that you should carefully study chapter 2, "Safety", and follow the instructions in this book.

OPERATING INSTRUCTIONS

System	POLYTRON®		
Type	PT 10-35 GT		
Issue	3 / 01.07.2018		
Page	4	of	32

1.1.1 RANGE OF VALIDITY

The information in these operating instructions relates to the POLYTRON® identified as follows:

Manufacturer: KINEMATICA AG, CH-6014 Luzern
 Brand name: POLYTRON®
 Product name: POLYTRON® PT 10-35 GT

Order No.	Identification
11010103	POLYTRON® PT 10-35 GT drive, 115V, with US-plug & extension for stand ST-P10/600
11010104	POLYTRON® PT 10-35 GT drive, 230V, with EU-plug & extension for stand ST-P10/600
11010105	POLYTRON® PT 10-35 GT drive, 230V, with CH-plug & extension for stand ST-P10/600
11010106	POLYTRON® PT 10-35 GT drive, 230V, with GB-plug & extension for stand ST-P10/600
11095000	adapter for dispersing aggregates of type PTA
11040050	stand ST-P10/600
11095002	vessel holder for ST-P10/600
35040006	stand ST-P20/600
11095001	extension Ø13 x 200 mm for stand ST-P20/600

1.1.2 TARGET AUDIENCE

These operating instructions are intended for all authorised users of our machines/equipment. We distinguish different user roles, taking account of the different demands placed on the user by the activity to be carried out.

You will find the definitions of user roles with the demands on the user in chapter 2, "Safety". You can fulfil one or more of these roles, provided that you meet the corresponding demands.

1.2 ORGANISATIONAL MATTERS

If you are unable to find the answer to any question in the operating instructions, please contact the equipment manufacturer directly.

1.2.1 LOCATION OF THE OPERATING INSTRUCTIONS

The operating instructions can only be of use to you if you always have them at hand. They should therefore always be kept at the place where the equipment is used.

1.2.2 MANUFACTURER CONTACT ADDRESS

KINEMATICA AG

Luzernerstrasse 147a
CH-6014 Lucerne

TEL: +41 41 259 65 65
FAX: +41 41 259 65 75
e-mail: laboratory@kinematica.ch

System	POLYTRON®		
Type	PT 10-35 GT		
Issue	3 / 01.07.2018		
Page	6	of	32

1.3 WARNING NOTICES

Please be aware of the meaning of the following warning signs:



SAFETY INSTRUCTIONS MUST BE OBSERVED TO ENSURE SAFE OPERATION .



THIS SYMBOL INDICATES HIGH VOLTAGE, WITH RISK TO HEALTH AND ENVIRONMENT.



**CAUTION!
BEWARE OF HOT SURFACE.**



**CAUTION!
DEVICE NOT DESIGNED FOR USE IN
EXPLOSION DANGER ENVIRONMENT.**

OPERATING INSTRUCTIONS			
System	POLYTRON®		
Type	PT 10-35 GT		
Issue	3 / 01.07.2018		
Page	7	of	32

2 SAFETY

This chapter is directed at all users of KINEMATICA laboratory equipment. It includes information on safe and optimum use.

2.1 SUMMARY

Any incorrect use of the installed equipment can be dangerous. Inadequately trained users can cause material damage and personal injury. This chapter informs you about the safety concept and the requirements for safe and optimum use of the equipment.

All those authorised to operate, service and repair the equipment are required to study chapter 2, "Safety".

2.2 SAFETY CONCEPT

The safety concept sets down the entitlement to use the equipment and the responsibilities of the individual users.

The machines and equipment are designed and constructed according to the state of the art and the recognised safety rules.

2.2.1 INTENDED USE OF THE EQUIPMENT

The equipment is designed and constructed for the following use:

- Dispersion and homogenisation of pumpable fluid products in accordance with the technical specifications (see point 3.5) and compatibility with the materials coming into contact with the products.

If you use the equipment for any purpose other than those listed, the manufacturer cannot be held liable for any resulting damage.

2.2.2 IMPROPER USE

Any use other than the "intended use" without the written approval of the manufacturer or any operation outside the technical limits of use is improper use.

2.2.3 USER ROLES

OPERATING INSTRUCTIONS

System	POLYTRON®		
Type	PT 10-35 GT		
Issue	3 / 01.07.2018		
Page	8	of	32

To guarantee safety, we place requirements on the users of the equipment that must be met without fail. Only persons meeting the requirements are authorised to work with the equipment.

We describe all those who work with the equipment as users. Since the requirements of these users are very much dependent on their activity, we distinguish the following user roles.

Contract partner:

The manufacturer can impose legal obligations on the contract partner when the equipment is purchased. The contract partner is obliged to ensure that the equipment is properly used.

Operating company:

The operating company ensures that the equipment is properly used and authorises persons who are entitled to work with the equipment in any one of the defined user roles. They are under the obligation to instruct the users.

Note:

Contract partner and operating company can be the same person.

Service technician:

The service technician is an employee of the operating company and looks after the equipment in special operating mode(s). He is a specialist with mechanical, electrical and electronic professional training. The service technician undertakes commissioning, decommissioning service and repair of the equipment. He must be appropriately trained to be able to carry out the service work required.

Operator:

The operator turns the equipment on and off. In the event of an alarm signal he informs the service technician.

OPERATING INSTRUCTIONS			
System	POLYTRON®		
Type	PT 10-35 GT		
Issue	3 / 01.07.2018		
Page	9	of	32

2.2.4 DANGER AREA

System/equipment

The system danger area includes the whole system/equipment including the connecting lead and controls.

Proximity danger area

This refers to all areas within a defined distance of the equipment.

User danger area

This danger area includes all persons working with the equipment.

2.2.5 AREAS OF RESPONSIBILITY

In order that the system/equipment can be used safely and without risk, the users in various roles bear the responsibility for particular danger areas.

Contract partner:

The contract partner bears the responsibility for the "proximity danger area".

Operating company:

The operating company bears the responsibility for the "user danger area". Only those users may be authorised to operate the system/equipment who fulfil all requirements of the user roles concerned. In doing so, attention must be paid to the following points:

- It is to be ensured that all users of the system/equipment have fully read and understood chapter 2, "Safety" and act accordingly in a safety-conscious manner.
- It is to be ensured that no unauthorised person carries out work with the system/equipment.
- It is to be ensured that users are informed of the possible risks and dangers connected with the system/equipment.
- It is to be ensured that those being trained or engaged in general training are under the permanent supervision of a trained and authorised person.

OPERATING INSTRUCTIONS

System	POLYTRON®		
Type	PT 10-35 GT		
Issue	3 / 01.07.2018		
Page	10	of	32

Service technician:

The service technician bears the responsibility for the "system/equipment danger area". He ensures that the system/equipment is at all times free from technical faults, safe and functions correctly.

2.2.6 GENERAL SAFETY RULES

Observe the following general safety rules:

- follow these operating instructions,
- in addition, observe the legal obligations and requirements for accident prevention and environmental protection of the country in which you operate the equipment,
- do not make any modifications to the equipment without the written authorisation of the manufacturer,
- only original replacement parts may be used for repairs,
- before any service work on the equipment, it must be ensured that the electrical supply is switched off,
- after any service, maintenance or repair work has been carried out on the system/equipment, it must be given a test run by the service technician.
- depending on the place at which it is installed, circumstances may require that hearing protection is worn when remaining in the vicinity of the equipment for long periods.

2.3 RESIDUAL DANGERS

When the system/equipment is used in accordance with rules and regulations, residual dangers are minimal.

Residual danger	Countermeasures
Tripping over feed or return lines	These should be laid appropriately.
Breakage of glass containers	Wear protective clothing (goggles etc.).  
Spitting of the product	
Hearing loss due to loud noise.	According to the application ear protection must be used. 
Tilting of the device	Use stable, non-slip base and the safety stands from KINEMATICA

System	POLYTRON®		
Type	PT 10-35 GT		
Issue	3 / 01.07.2018		
Page	11	of	32



IN EVERY CASE THE ELECTRICAL INSTALLATION HAS TO BE DONE BY TECHNICIAN!

2.4 WARNINGS

	• Ensure that the rated voltage of the equipment matches the supply. • Before changing any dispersing aggregate, the line cord has to be plugged out • When a electrical power blackout occurs, the device has to be switched- off using the main switch, avoiding the device to restart automatically. • IT IS IMPORTANT THAT THE MAINS SUPPLY WHERE THE DEVICE IS PLUGGED IN COMPLIES WITH THE INFORMATION ON THE TYPE LABEL AND THE INTERNATIONAL STANDARDS FOR POWER SUPPLIES. IF NOT, SUCCESSFUL OPERATION CANNOT BE GUARANTEED
	• In the event that hazardous chemicals or materials that endanger health can influence the surroundings or use of the equipment, appropriate countermeasures must be taken.
	• At long term use the aggregate and the coupling may get hot – danger of skin burn.
	• The equipment may not be operated in explosive areas • It is not allowed to work with fluids which are highly inflammable. • It is not allowed to mix materials which can cause strong exothermal reactions

WARNINGS: continued

OPERATING INSTRUCTIONS

System	POLYTRON®		
Type	PT 10-35 GT		
Issue	3 / 01.07.2018		
Page	12	of	32

	• THE DEVICE MAY ONLY BE OPENED BY AUTHORISED KINEMATICA SERVICE STATIONS. • Ensure that enough free space is available at the backside of device, so that effective air flow and cooling is assured. Insufficient cooling may lead to a decrease of power output. • The device has to be placed in a manner that dirt or fluids cannot penetrate through the ventilation slots at back side of the drive. • POLYTRON® dispersion aggregates may not be operated dry – the lower sleeve bearing is cooled and lubricated by the medium being processed. Running dry will destroy the sleeve bearing. • The dispersing aggregates should be cleaned after every operation. • Never pull the coupling during operation – the aggregate could fall out of the coupling.
	• When the line cord is plugged, never touch the saw teeth of the aggregate – danger of injuries due to rotating shafts and blades
	KINEMATICA AG products comply with all the usual CE directives, carry the CE marking and are delivered with a corresponding declaration of conformity.

System	POLYTRON®		
Type	PT 10-35 GT		
Issue	3 / 01.07.2018		
Page	13	of	32

3 DESCRIPTION OF THE EQUIPMENT

The PT 10-35 GT system conforms to all current and relevant CE regulations, carries the CE sign and is supplied with the relevant certificate of conformity.

3.1 OVERVIEW

The POLYTRON® System PT 10-35 GT prepares the finest dispersions and homogenisations and is suitable for use with working volumes up to 5000 ml (depending on the viscosity of the product and the dispersing aggregate used). The drive can be equipped with a wide range of POLYTRON® dispersing-aggregates. The aggregates are easy to change using the quick-coupling.

Using the adapter (optional, order no.: 11095000), PTA-aggregates can also be equipped.


For a ready-to-use system you need:

- Drive PT 10-35 GT
- A stand, ST-P10/600 or ST-P20/600
- A dispersing aggregate
- A mains connection according to the type label

On the following pages pictures and a description of the main components of the complete system PT 10-35 GT can be found.

OPERATING INSTRUCTIONS

System	POLYTRON®		
Type	PT 10-35 GT		
Issue	3 / 01.07.2018		
Page	14	of	32

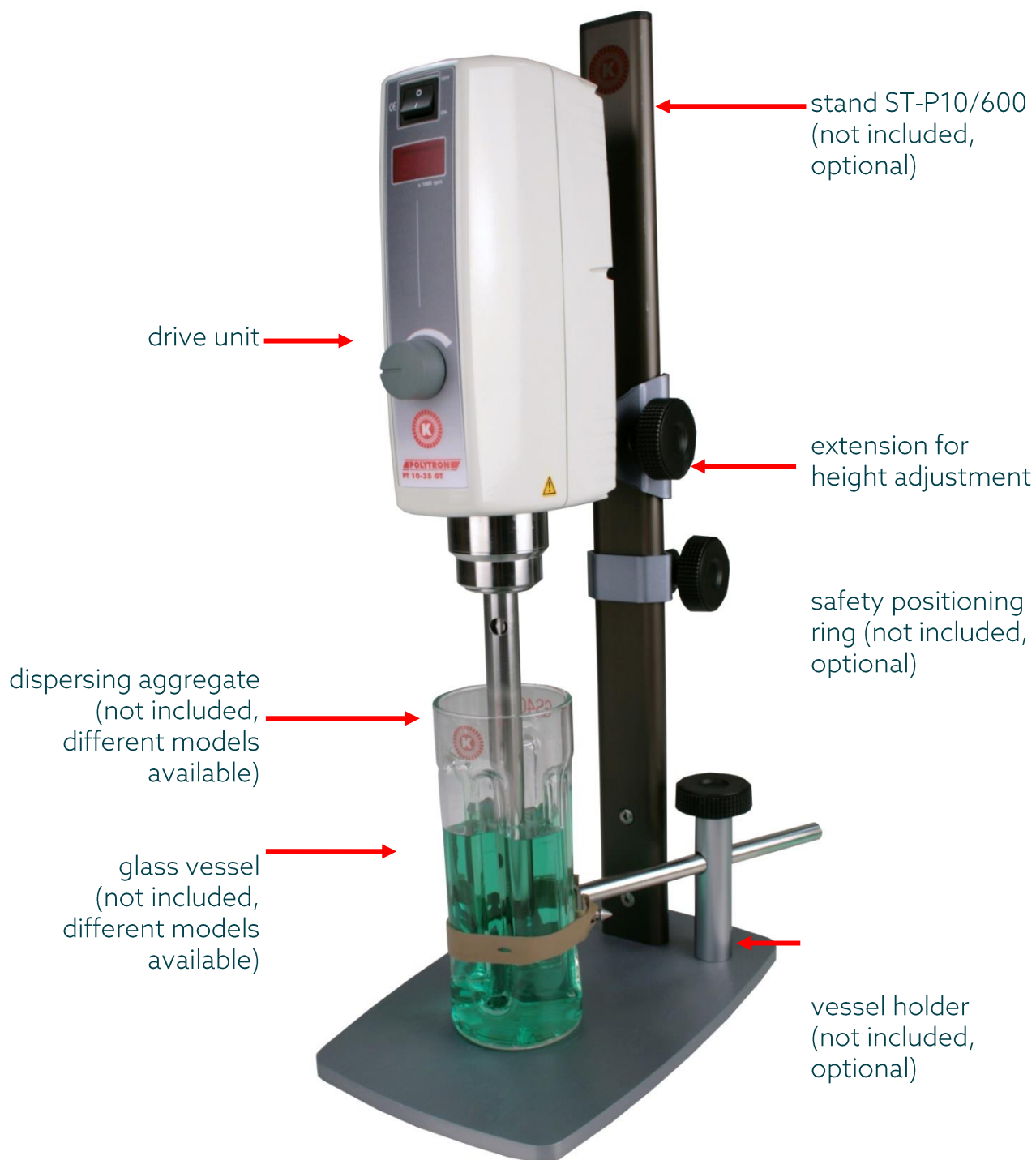

3.2 DRIVE UNIT

The drive unit PT 10-35 GT is equipped with a 1'200 W motor and will be started using the main switch placed at the front side. The speed will be adjusted using the turning


OPERATING INSTRUCTIONS

System	POLYTRON®		
Type	PT 10-35 GT		
Issue	3 / 01.07.2018		
Page	15	of	32


knob on the front side of the device. Within the speed range of up to 30'000 rpm, the integrated closed loop speed control will tune the speed to any adjusted value, independent from load changes. The actual speed can be read from the digital display (LED)

Furthermore the electronic control system is equipped with safety functions with automatic shut down to prevent the drive being overheated or blocked. A detailed description of possible errors and countermeasurements is listed in chapter 6 „TROUBLE SHOOTING“

BACK SIDE

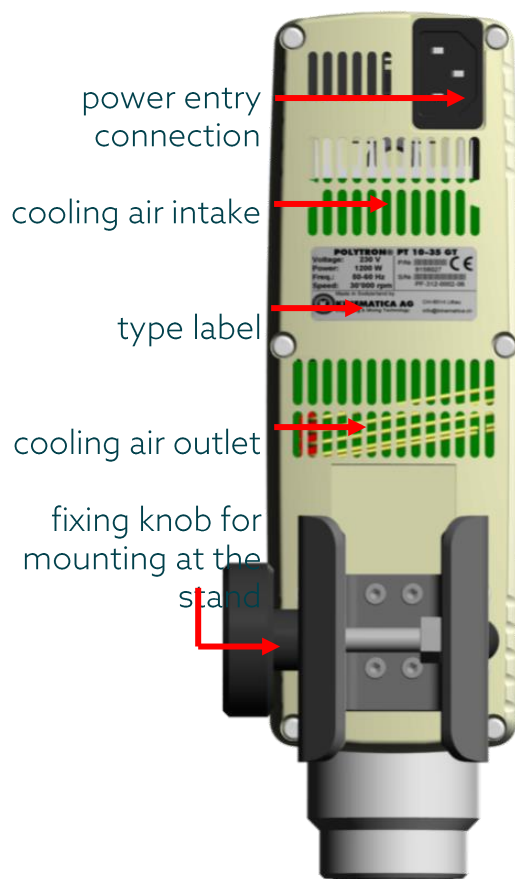

FRONT SIDE

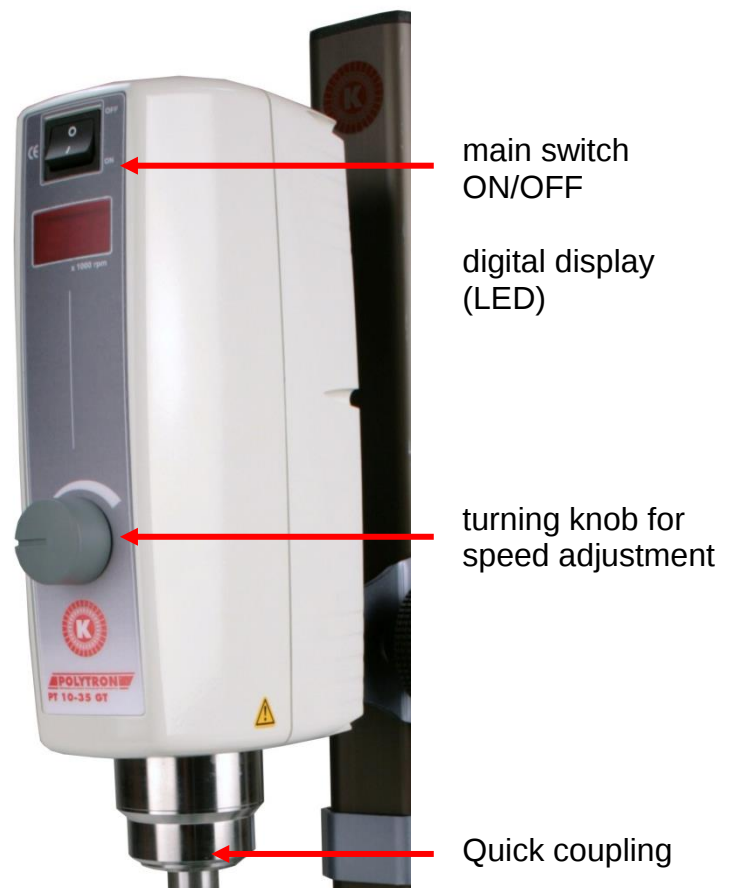


System	POLYTRON®		
Type	PT 10-35 GT		
Issue	3 / 01.07.2018		
Page	16	of	32

3.3 STANDS ST-P10/600 & ST-P20/600

The drive PT 10-35 GT is designed for using with a stand. KINEMATICA recommends the safety stands ST-P10/600 or optional ST-P20/600. According to the size of the working vessel a vessel holder may be used.

Description of the main components of the two stands:

stand ST-P10/600

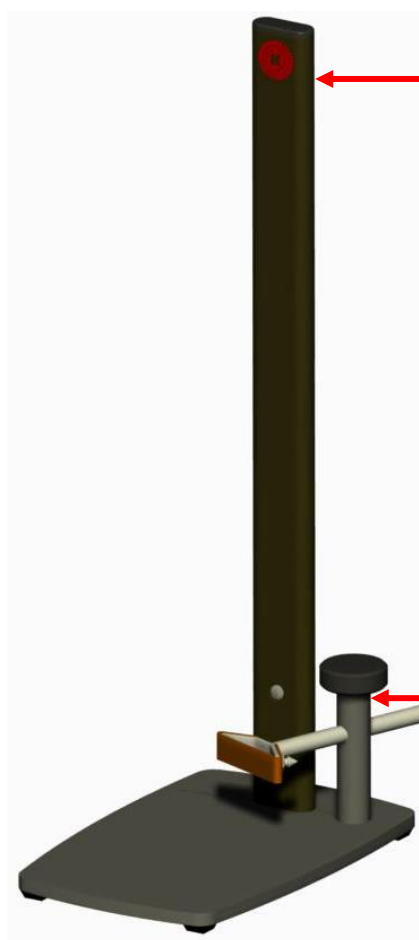

column with scratch-resistant surface made of aluminium


column made of stainless steel

boss head for extension Ø13 mm

vessel holder

anodised base plate

stand ST-P20/600

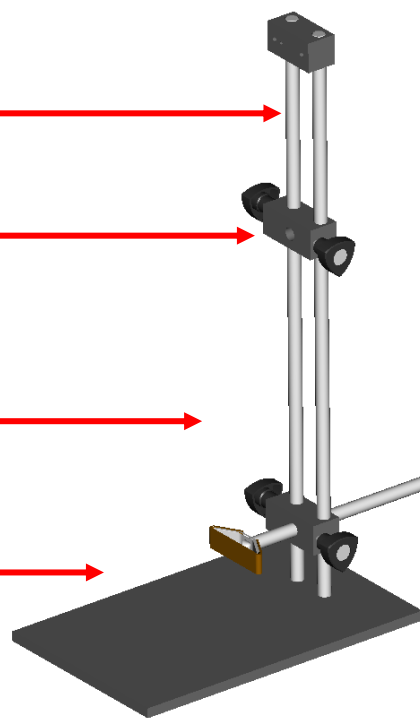

vessel holder

System	POLYTRON®		
Type	PT 10-35 GT		
Issue	3 / 01.07.2018		
Page	17	of	32

3.4 POLYTRON® DISPERSING-AGGREGATES


A wide range of dispersing aggregates with diameters of 5, 7, 12, 20, 30, 36 and 40 mm are available. For questions concerning POLYTRON® dispersing aggregates please get in contact with KINEMATICA or your local specialist dealer.

Due to the quick coupling, POLYTRON® dispersing aggregates can be fitted and replaced very easily. The three pictures below show the mounting of an aggregate.

For easy mounting, the crown head of the aggregate and the coupling should be free of dirt.

Push the slider of the coupling upwards, at the same time push aggregate into the coupling.

Push the aggregate upwards and at same time twist it a little until it has snapped in and the slider returns downwards

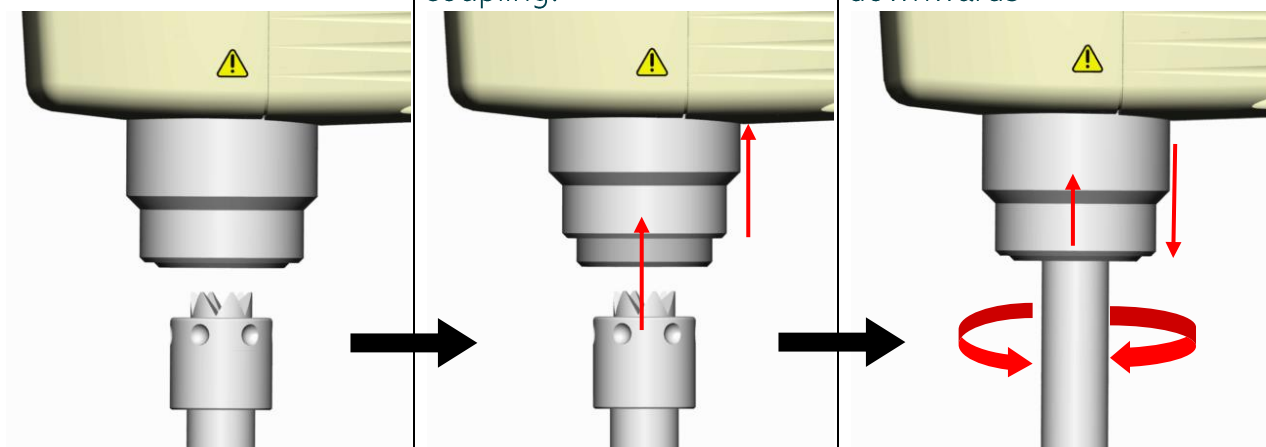

To remove the aggregate, proceed in reverse direction.




CAUTION !:

BEFORE PUSHING UP THE SLIDER OF THE COUPLING, HOLD YOUR DISPERSING AGGREGATE IN ONE HAND AVOIDING THE AGGREGATE TO FALL DOWN WHEN PUSHING UP THE SLIDER.

System	POLYTRON®		
Type	PT 10-35 GT		
Issue	3 / 01.07.2018		
Page	18	of	32



CAUTION!
WHEN REPLACING AN AGGREGATE
THE LINE CORD HAS TO BE
DISCONNECTED FROM THE POWER
SUPPLY

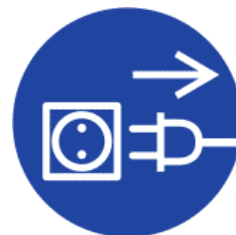

CAUTION!
AT LONGTERM USE AT HIGH LOADS, THE AGGREGATE
& THE COUPLING CAN HEAT UP TO HIGH
TEMPERATURE LEVELS - DANGER OF SKIN BURN .

3.5 TECHNICAL SPECIFICATIONS

Drive PT 10-35 GT	
motor type	universal motor
supply voltage	3.5.1.1.1.1.1 2 100...120 V ~
supply frequency	50/60 Hz
max. speed	3.5.1.1.1.1.2 30'000 rpm
Speed indication	3.5.1.1.1.1.3 LED, 3 digits, actual speed
Speed measurement	3.5.1.1.1.1.4 Digital, direct by HAL-probe (no calibration needed)
input power	3.5.1.1.1.1.5 Max. 1'200 W
output power	3.5.1.1.1.1.6 Max. 800 W
soft-start	Yes
noise emission (drive only)	lower than 75 dB(A) at 30,000 rpm
direction of rotation	clockwise, seen from above
ambient temperature	0 - 40°C
relative humidity	95% max.
certified standards	EMC according to: IEC/EN 61000-6-3 safety according to: IEC/EN 61000-6-2 IEC/EN 61010-2-51
protection type	IP 20
max. period of continuous operation	depending on load & speed 100 % duty time is possible.
Dimensions	see drawing
weight	drive unit 3.2 kg

OPERATING INSTRUCTIONS

System	POLYTRON®		
Type	PT 10-35 GT		
Issue	3 / 01.07.2018		
Page	19	of	32

Stand PT-ST P10/600

dimensions HxBxT	600x200x300 mm
weight	ca. 2.5 kg

Stand PT-ST P20/600

dimensions HxBxT	600x200x315 mm
weight	ca. 3.8 kg

4 INSTALLATION

4.1 UNPACK

Open the dispatch box and check that the contents agrees with the delivery note.




CHECK ALL PARTS FOR POSSIBLE TRANSPORT DAMAGE. INFORM US OR YOUR DEALER IMMEDIATELY ABOUT ANY DISAGREEMENT OR FAULT. If possible SEND US DIGITAL PHOTOS BY EMAIL TO LABORATORY@KINEMATICA.CH

System	POLYTRON®		
Type	PT 10-35 GT		
Issue	3 / 01.07.2018		
Page	20	of	32

4.2 ASSEMBLING OF THE DRIVE


The PT 10-35 GT drive arrives fully equipped with the extension suitable for the stand ST-P10/600. Just mount the drive at the stand and plug the line cord into the power inlet connection at the back side. When the double rod stand ST-P20/600 will be used, a suitable extension can be attached to the drive. The de/assembling of these components is explained below.

Dis/assembling of the extension for the stand ST-P10/600

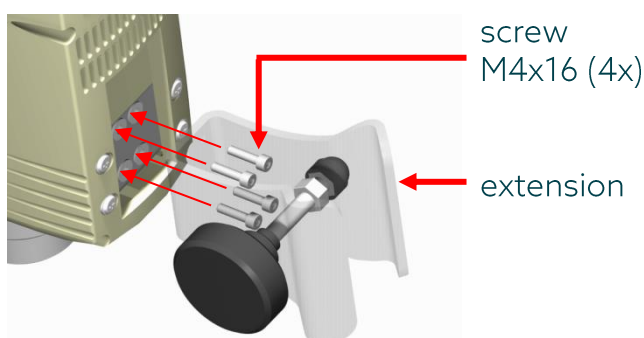

For de/assembling, please use the supplied tool.

Dis/assembling of the extension Ø13x200 (optional) for the stand ST-P20/600

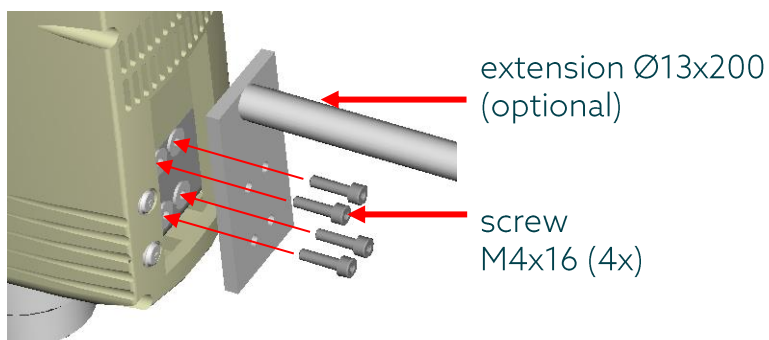

For de/assembling, please use the supplied tool.




CAUTION!
DO NOT USE OTHER SCREWS. SCREWS WHICH ARE TOO LONG MAY DAMAGE INTERNAL COMPONENTS OF THE DRIVE.

4.3 ASSEMBLING OF THE STAND WITH ACCESSORIES

Assembling of the stand ST-P10/600


OPERATING INSTRUCTIONS

System	POLYTRON®		
Type	PT 10-35 GT		
Issue	3 / 01.07.2018		
Page	21	of	32

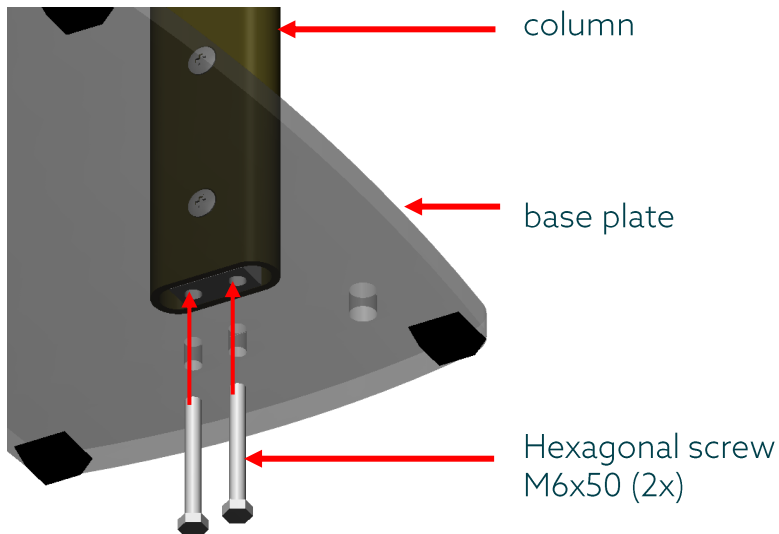


Attach the column to the base plate using the two screws. Fix the screw using the provided tool

Assembling of the vessel holder for ST-P10/600

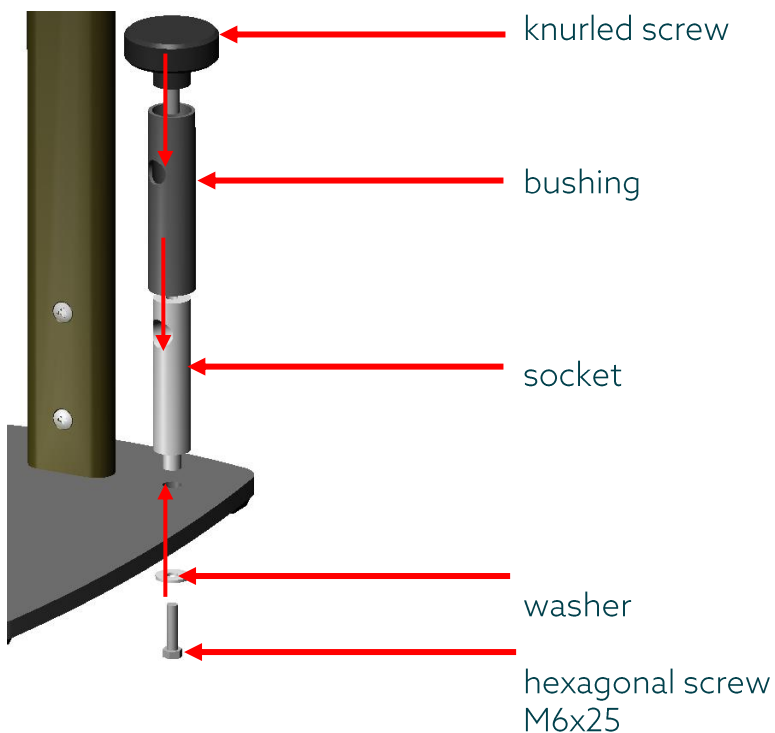

Insert the vessel holder extension as shown on the picture below

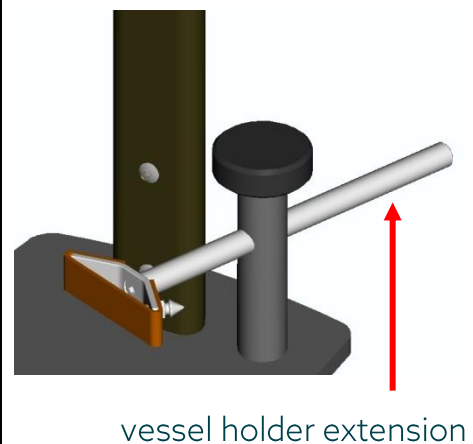

4.4 STARTING UP

- Ensure that the voltage marked on the type label complies with your power supply

OPERATING INSTRUCTIONS

System	POLYTRON®		
Type	PT 10-35 GT		
Issue	3 / 01.07.2018		
Page	22	of	32

- Immerse the aggregate in the product. The optimum depth of immersion of the dispersion attachment is about 70 % of the total depth of product in the container. Never immerse the upper coupling head and upper hole of the dispersion aggregate.
- Turn on at the main switch.
- Set the desired speed using the turning knob

Remark:

In some cases it can be advantageous to place the vessel excentric in order to get more turbulences for better mixing of the product.



POLYTRON® DISPERSING AGGREGATES MAY NOT BE OPERATED DRY - THE LOWER SLEEVE BEARING IS COOLED AND LUBRICATED BY THE MEDIUM BEING PROCESSED. RUNNING DRY WILL DESTROY THE SLEEVE BEARING.

THE DISPERSING AGGREGATES SHOULD BE CLEANED AFTER EVERY OPERATION.

NEVER TOUCH THE ROTATING AGGREGATE - DANGER OF SEVERE CUT-INJURIES.

System	POLYTRON®		
Type	PT 10-35 GT		
Issue	3 / 01.07.2018		
Page	23	of	32


5 MAINTENANCE

Your POLYTRON® unit is designed for ease servicing. Nevertheless, it is essential to inspect your equipment carefully and to clean it thoroughly. Drawings of the separate components are to be found in the appendix.



THE EQUIPMENT MUST BE DIS-CONNECTED FROM THE ELECTRICAL SUPPLY:

- DURING ANY WORK ON THE EQUIPMENT, IN ORDER TO AVOID ANY PERSONAL INJURY OR OTHER DAMAGE
- WHEN CHANGING OR REMOVING THE DISPERSING AGGREGATE

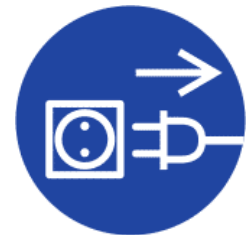

5.1 DRIVE

Under normal working conditions, the PT 10-35 GT drive requires no servicing. Parts such as the motor brushes and bearings are subject to natural wear.

The following are standard values for wear parts of the drive:

- bearing: several 1000's of operating hours.
- brushes: Approx. 150 hours of operating at rated power.

We strongly recommend that service work and repairs should be carried out only by authorised KINEMATICA service centres or by KINEMATICA directly, where original replacement parts are available.

Any unauthorised modification or manipulation of the unit or its equipment leads to immediate annulment of the warranty.

System	POLYTRON®		
Type	PT 10-35 GT		
Issue	3 / 01.07.2018		
Page	24	of	32

5.2 CLEANING AND MAINTENANCE OF THE AGGREGATES


EC-aggregates, (EC stands for EASY CLEAN) are easy to disassemble and clean. With some restrictions(*) they can be sterilised in an autoclave by several methods. In the following three chapters the de/assembling procedures will be described.

(*) Depending on the frequency and the intensity of the sterilisation process, the life span of the bearings may be reduced due to loss of grease during evacuation in a autoclave.



CAUTION!
BE AWARE OF SHARP EDGES AND TEETH WHEN DE/ASSEMBLING AGGREGATES – DANGER OF CUT-INJURIES


5.2.1 Dis/assembling of EC-aggregates with size Ø5 and Ø7 mm

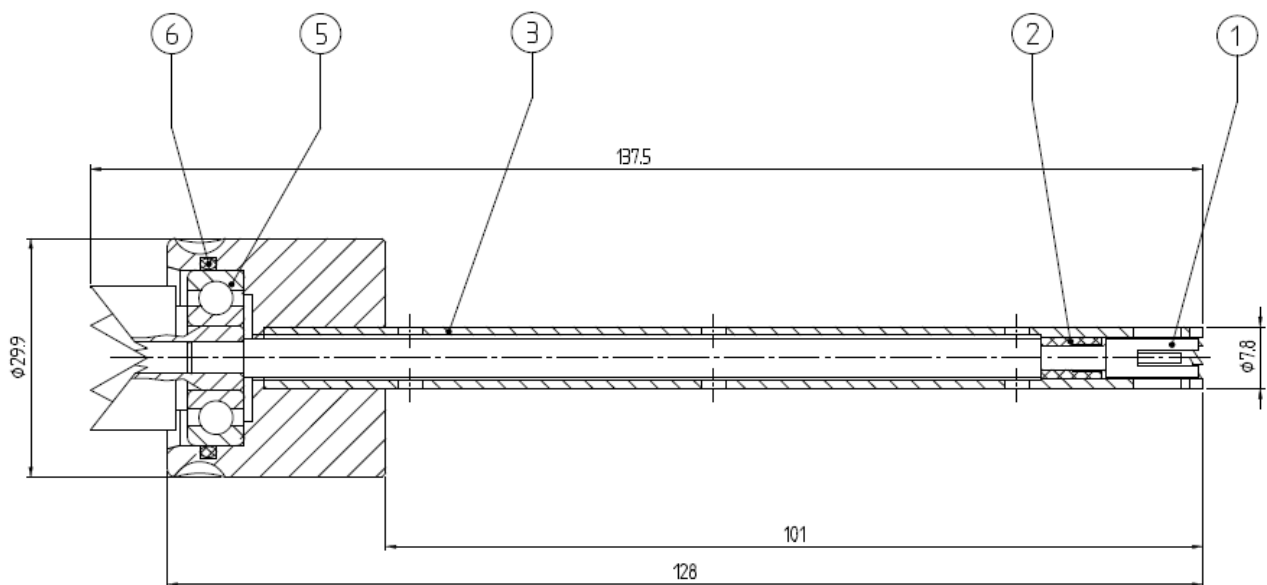

1. Push the shaft (1) from the right side out of the stator pipe.
2. Remove the slide bearing(2) from the shaft.
3. Remove the o-ring(6).
4. After the cleaning process, reassemble the aggregate in reverse order.


System	POLYTRON®		
Type	PT 10-35 GT		
Issue	3 / 01.07.2018		
Page	25	of	32

5.2.2 Dis/assembling of EC-aggregates with size Ø 12 und Ø 20 mm

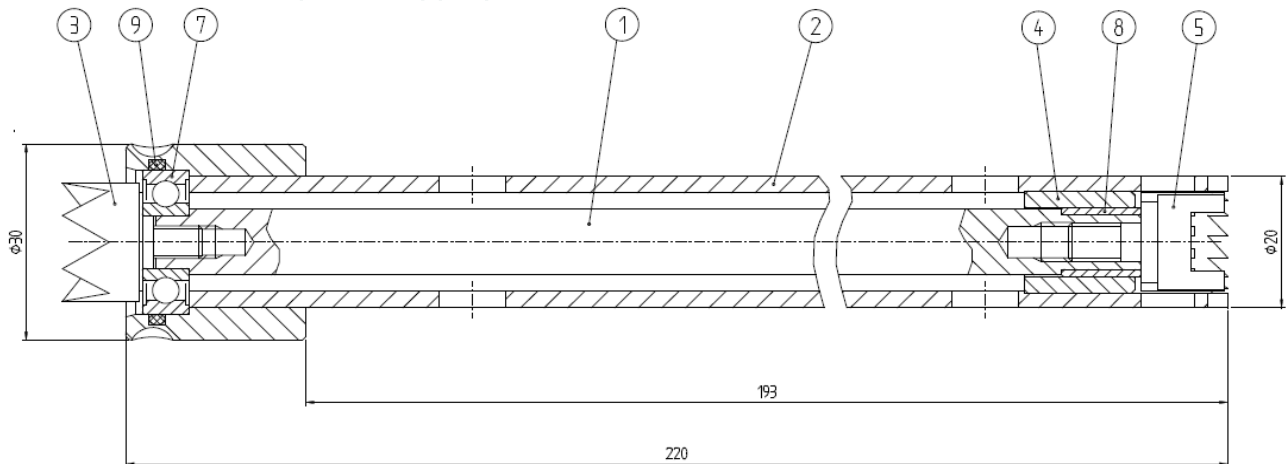

1. Fix the crown head(3) using the tool. Unscrew the rotor (5) using a suitable tool. Pull out the shaft(1)of the stator tube.
2. Remove the o-ring(9) out of the groove.
3. Unscrew the crown head from the shaft.
4. After the cleaning process, reassemble the aggregate in reverse order.

5.2.3 Dis/assembling of EC-aggregates with size Ø30+

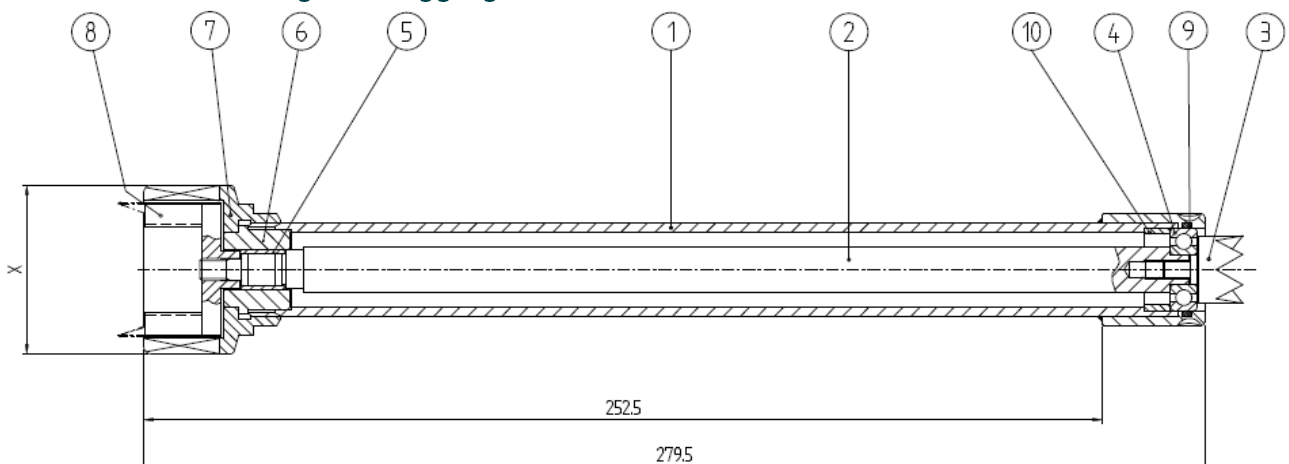

1. Fix the crown head (3) using the tool. Unscrew the rotor (8) using a suitable tool. Pull out the shaft (2) of the stator tube.
2. Unscrew the stator(7).
3. Pull out the shaft (2) of the stator tube(1).
4. Remove the o-ring (9) out of the groove.
5. Unscrew the crown head(3).
6. After the cleaning process, reassemble the aggregate in reverse order.

OPERATING INSTRUCTIONS

System	POLYTRON®		
Type	PT 10-35 GT		
Issue	3 / 01.07.2018		
Page	26	of	32

5.2.4 Criteria for replacement of the ball bearings

Basically, it is the operating company that determines when and how often the bearings are to be changed. They should, however, be changed at the latest if:

- rotor and stator are touching
- there is an increase in vibration,
- the rotor projects above the edge of the stator.

For a replacement of the ball bearings please contact a certified KINEMATICA service center or KINEMATICA directly.

Any unauthorised modification or manipulation of the unit or its equipment leads to immediate annulment of the warranty.

6 TROUBLE SHOOTING

PROBLEM	REASON	CORRECTIVE MEASURES
Unusual noises	Damaged drive bearings	Change ball bearings. Trace & replace defective parts (shaft, bearings)
	Damaged aggregate bearings	Change bearings
	Rotor/stator interference	Trace and replace defective parts (shaft, bearings)
Drive automatically stops & an alternating beeping sound is audible and the message "HO" is shown on the display	Inadequate ventilation	Check if ventilation slots are clear. Cool down the device & restart
	Thermal overload	Check manner of use. Cool down the device & restart.
Vibrations	Bent shaft	Replace shaft
	Worn bearing(s)	Replace bearing(s)
	Defective coupling	Trace and replace defective parts
Drive does not start respectively suddenly stops. A constant beeping sound is audible and the message "BL" is shown on the display	Defective internal control	Send the equipment to the nearest authorised KINEMATICA service centre or directly to KINEMATICA AG
	Drive is blocked	Check the aggregate for solid particles which may be blocking the rotor, remove the particles, turn off the drive and restart operation.

OPERATING INSTRUCTIONS

System	POLYTRON®		
Type	PT 10-35 GT		
Issue	3 / 01.07.2018		
Page	27	of	32

7 ACCESSORIES

A large selection of special homogenising vessels is available, which can improve the efficiency of the processing of your product. Closed vessels to withstand vacuum or pressure can also be supplied.

Ask your dealer or contact KINEMATICA AG directly.

8 WARRANTY

KINEMATICA AG guarantees that the equipment that it produces will run free of any fault related to materials or manufacturing faults for **24 months**.

ADDITIONAL WARRANTY FOLLOWING REGISTRATION

Consumers who complete a warranty registration within 20 days of receiving their device will benefit from a free warranty extension of a further 12 months. This means that the total warranty period after registration is 36 months.

REGISTER NOW



kinematica.ch/warranty-upgrade

If thorough testing shows a fault to be due to either of the above causes, KINEMATICA AG guarantees that the equipment will be repaired or replaced free of charge. The guarantee does not cover parts that are subject to normal wear. It is void if any person other than an employee of KINEMATICA AG or their appointed representative has made modifications to the equipment or if the damage is due to failure to comply with the operating instructions, to carelessness, accident, incorrect use or incorrect supply voltage.

Please check our further warranty conditions and the correct procedure for a repair application on our website:

<http://www.kinematica.ch/en/contact-service/service-and-repair.html>

KINEMATICA AG reserves the right to make technical changes to the equipment without modifying equipment delivered earlier in the same way.

In the event of technical problems, for spare parts requirements or for advice, contact our regional appointed agent, your preferred dealer or us directly at:

KINEMATICA AG
Luzernerstr. 147a
CH-6014 Luzern
Switzerland

Tel. +41-41-259 65 65
Fax +41-41-259 65 75
Email service@kinematica.ch

OPERATING INSTRUCTIONS

System	POLYTRON®		
Type	PT 10-35 GT		
Issue	3 / 01.07.2018		
Page	29	of	32

9 DISPOSAL


The symbol of the crossed refuse container signifies that the product in the European Union must be supplied to separate collection. Labeled products must not be disposed with household waste, rather must be left at a collection point for recycling electrical and electronic equipment. Recycling helps to reduce the consumption of raw materials and to protect the environment.



System	POLYTRON®		
Type	PT 10-35 GT		
Issue	3 / 01.07.2018		
Page	30	of	32

Appendix A. dimensional drawings

Drive PT 10-35 GT

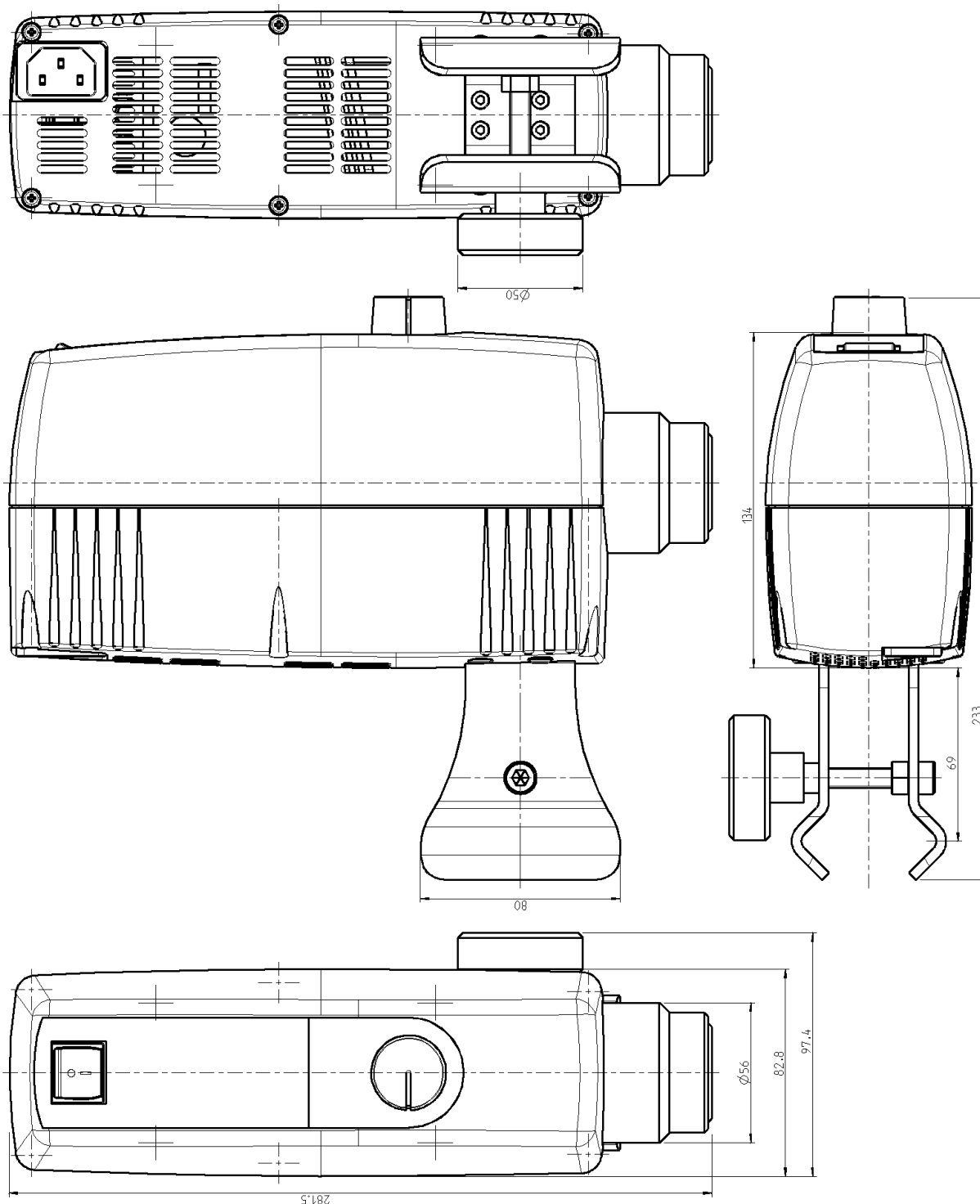

Stand ST-P10/600




KINEMATICA

Homogenizing perfected.

OPERATING INSTRUCTIONS

System	POLYTRON®		
Type	PT 10-35 GT		
Issue	3 / 01.07.2018		
Page	31	of	32

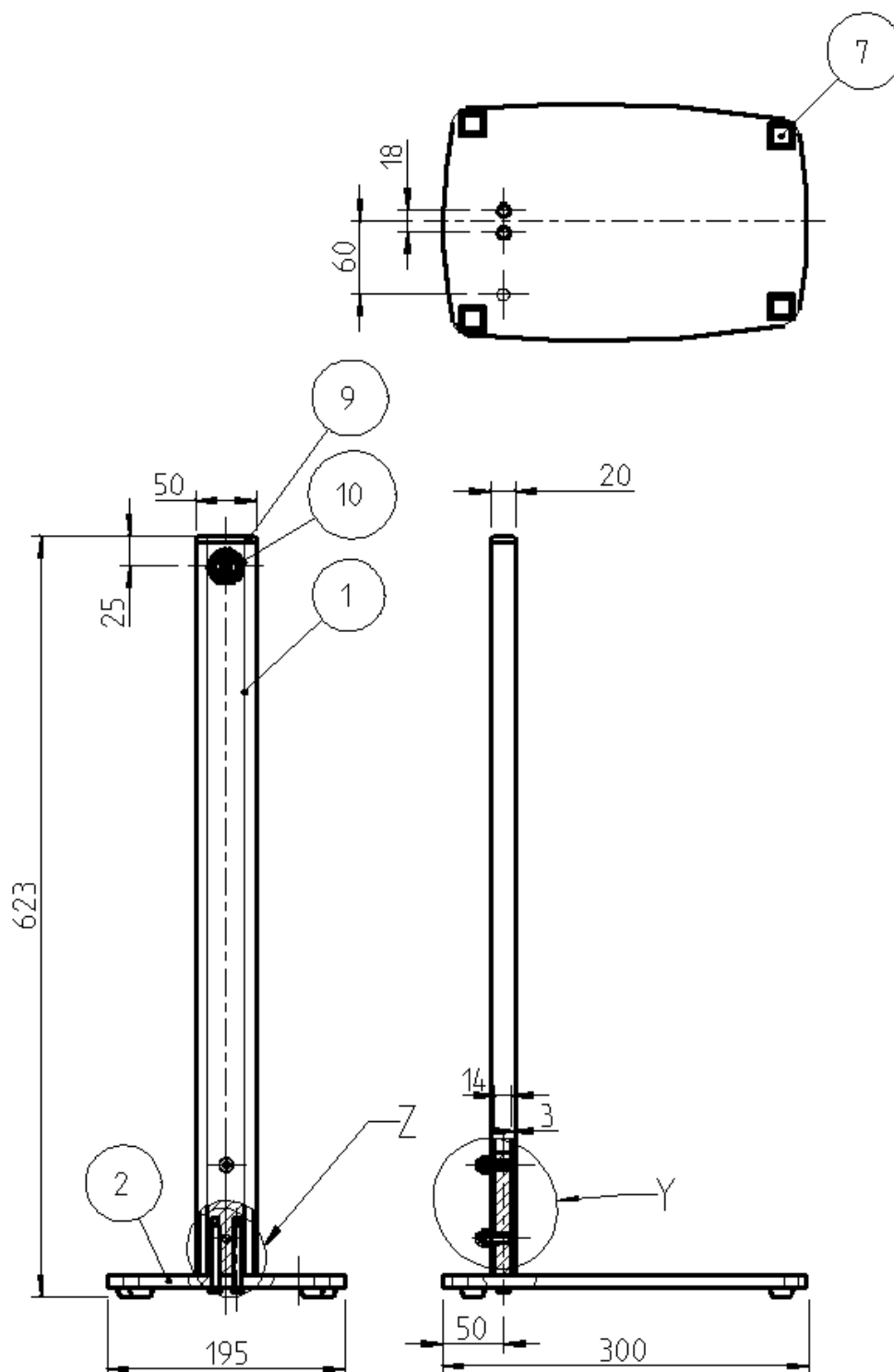

OPERATING INSTRUCTIONS

System	POLYTRON®		
Type	PT 10-35 GT		
Issue	3 / 01.07.2018		
Page	32	of	32

Appendix B. bill of materials

Assembly Group	Main Identification	Order No.	Production-No.	Product-Identification
Drives	Drive POLYTRON® PT 10-35 GT, 115 V	11010103	9158158	POLYTRON® PT 10-35 GT, 115 V
			9341277	Line cord with US-plug
	Drive POLYTRON® PT 10-35 GT, 230 V/EU	11010104	9158159	Drive POLYTRON® PT 10-35 GT, 230 V
			9341278	Line cord with EU-plug
	Drive POLYTRON® PT 10-35 GT, 230 V/CH	11010105	9158159	Drive POLYTRON® PT 10-35 GT, 230 V
			9341282	Line cord with CH-plug
Stands	Stand ST-P10/600	11040050	9158159	Drive POLYTRON® PT 10-35 GT, 230 V
			9341279	Line cord with UK-plug
	Stand ST-P20/600	35040006	9115323	ST-P10/600
			9341391	Flat spanner 10 mm
			9115268	ST-P20/600
			9900738	Cable holder
Options	Vessel holder for stand ST-P10/600	11095002	9900046	Screw M4x6
			9770007	Set of tools
			9754230	Vessel holder
	Adapter for PTA-aggregates	11095000	9135040	Vessel holder extension
			9800064	Rubber band
			9112088	Adapter PTA/PT-DA
	extension Ø13 x 200 mm for stand ST-P20/600	11095001	9754043	Extension Ø13x200
			9920416	Screw M4x16 (4x)
			9300112	Hex wrench